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
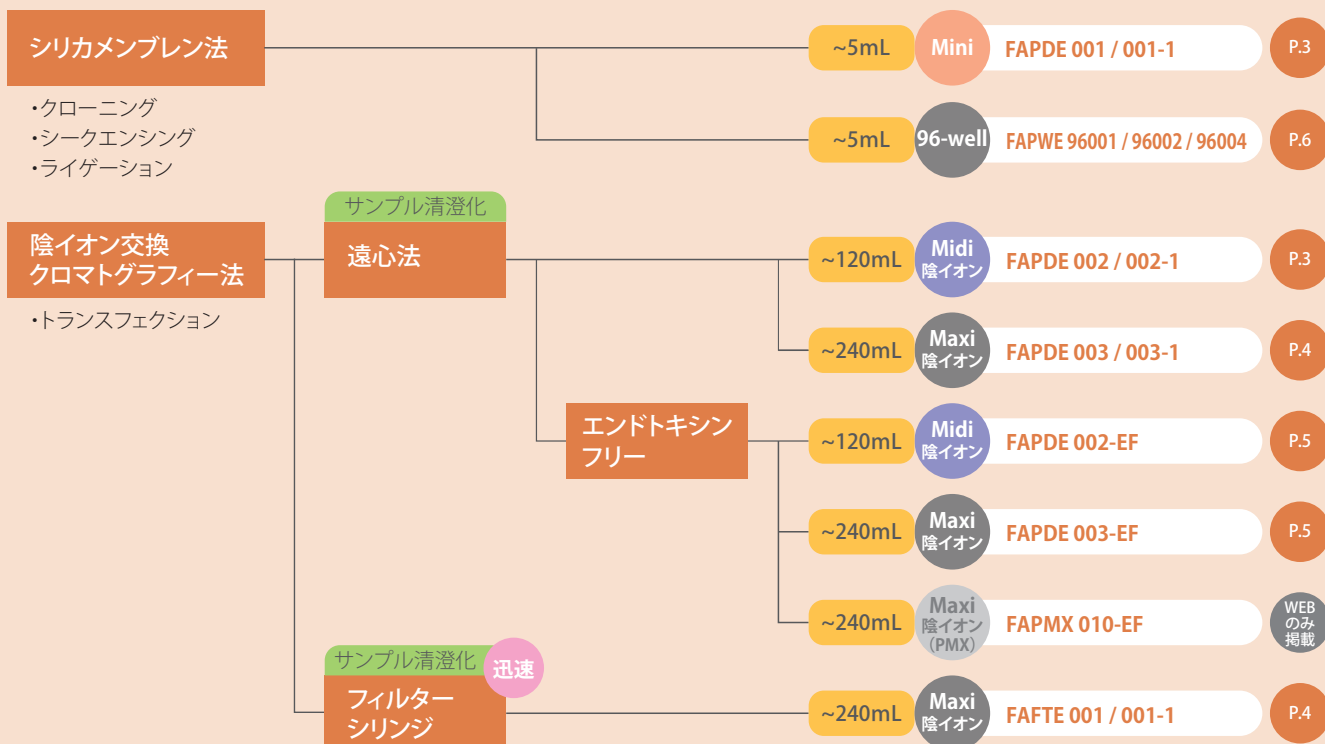
FAPD1 Buffer, FAPD2 Buffer, FAPD3 Buffer, WF Buffer (concentrated), Wash Buffer (concentrated), Elution Buffer, RNase A (liquid), FAPD Column, Collection Tube

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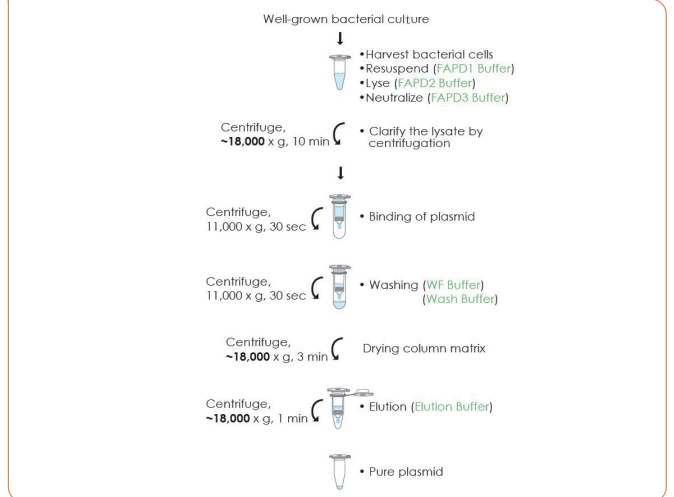FavorPrep Plasmid Extraction Midi Kit

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
PEQ Buffer, PM1 Buffer, PM2 Buffer, PM3 Buffer, PW Buffer, PEL Buffer, RNase A (liquid), PM Midi Column

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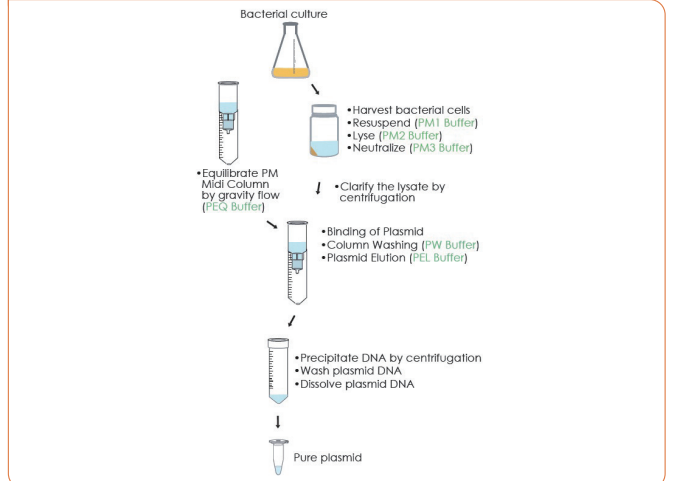
PEQ Buffer, PM1 Buffer, PM2 Buffer, PM3 Buffer, PW Buffer, PEL Buffer, RNase A (liquid), PM Maxi Column

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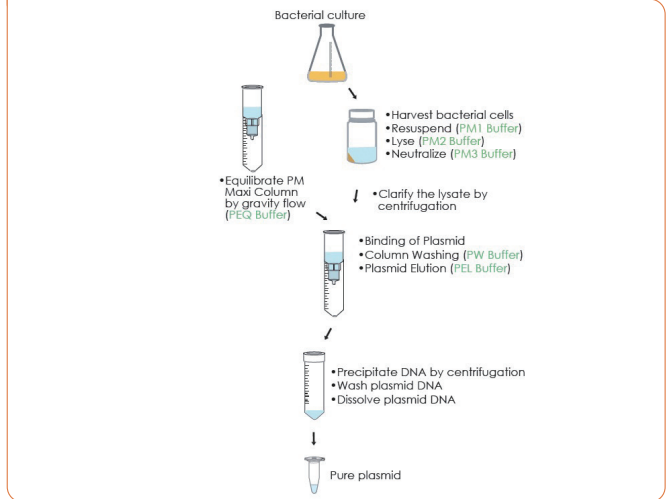
PEQ Buffer, PM1 Buffer, PM2 Buffer, PM3 Buffer, PW Buffer, PEL Buffer, RNase A (liquid), FavorFilter Maxi Cartridge, PM Maxi Column

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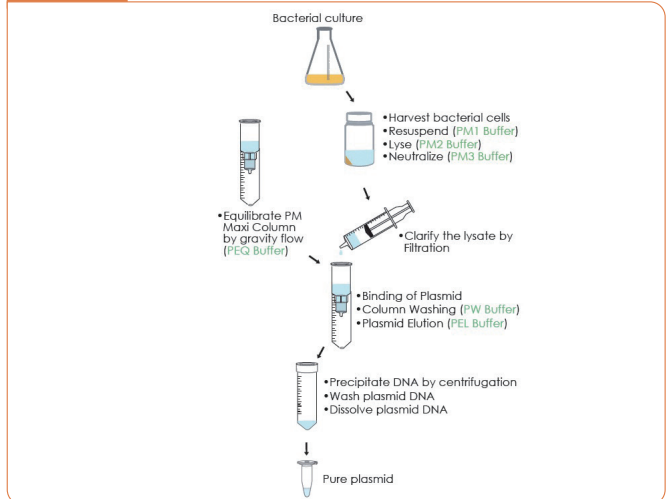FavorPrep Endotoxin Free Plasmid Extraction Midi Kit

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
PEQ Buffer, PM1 Buffer, PM2 Buffer, PM3 Buffer, PTR Buffer, PW Buffer, PEL Buffer, RNase A (liquid), PM Midi Column

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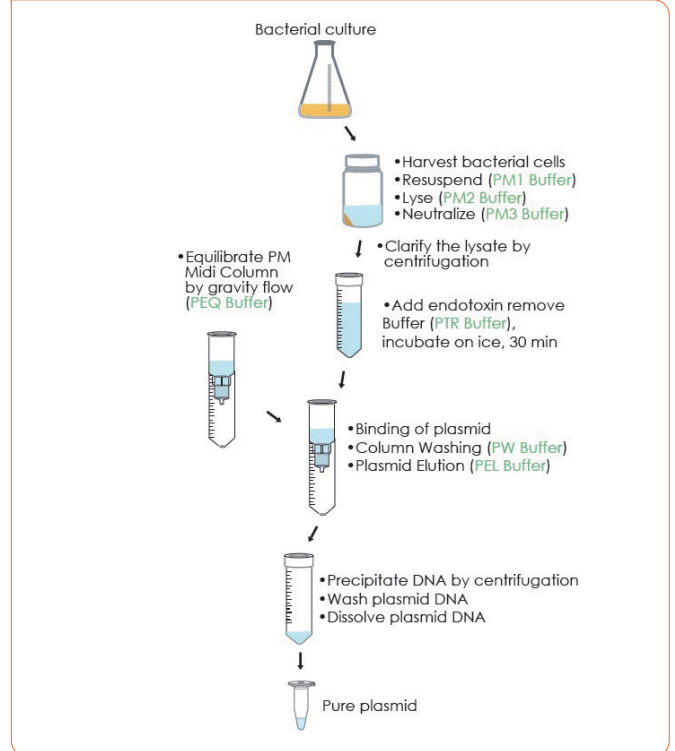FavorPrep Endotoxin Free Plasmid Extraction Maxi Kit

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
PEQ Buffer, PM1 Buffer, PM2 Buffer, PM3 Buffer, PTR Buffer, PW Buffer, PEL Buffer, RNase A (liquid), PM Maxi Column

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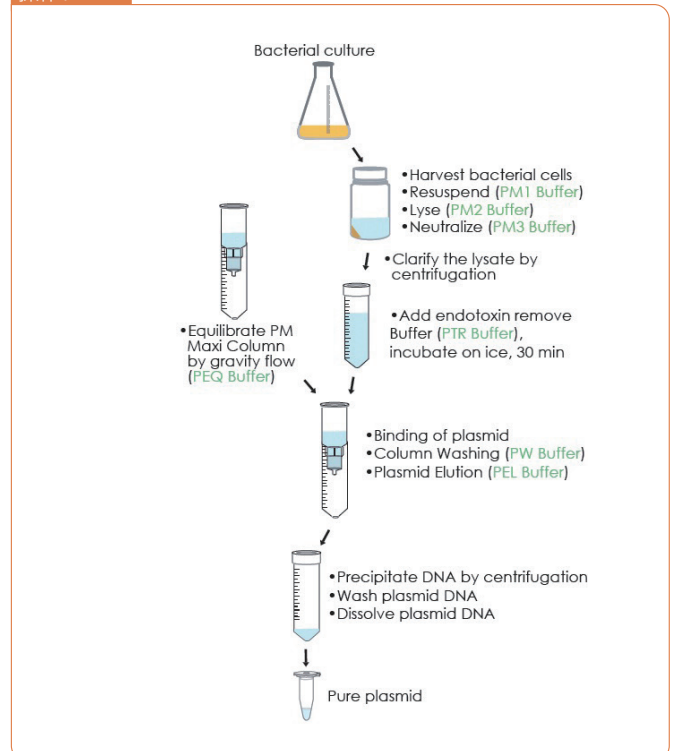FAPD1 Buffer, FAPD2 Buffer, FAPD3 Buffer, Wash Buffer (concentrated), Elution Buffer, RNase A (liquid), Filter Plate (96-Well Plasmid Binding plate), Collection Plate (96-Well 2 ml Plate), Elution Plate (96-Well PCR plate), Adhesive Film

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STEP 1. Collect bacterial cells and resuspend the cells

- Transfer bacterial cells to a Collection Plate (first Collection Plate)
- Seal with a Adhesive Film, and centrifuge at 5,600-6,000 x g for 3 min
- Add PDE1 Buffer
- Resuspend by pipetting

STEP 2. Lysis

- Add PDE2 Buffer
- Mix gently by pipetting 5 times

STEP 3. Neutralization

- Add PDE3 Buffer
- Mix by pipetting

STEP 4. Clarify lysate

- Seal with a Adhesive Film, and centrifuge at 5,600-6,000 x g for 10 min

STEP 5. Bind plasmid to Filter Plate:

vacuum processing

- Transfer the supernatant to Filter plate.
- Apply -12 inches Hg vacuum until the wells have emptied.

centrifuge processing

- Transfer the supernatant to Filter plate.
- Centrifuge at 4,500-6,000 x g for 2 min.

STEP 6. Wash the Filter Plate and dry the membrane of the Filter Plate

- Add Wash Buffer.
- Apply vacuum at -12 inches Hg for 2 min.
- Top the Filter Plate lips on paper towel
- Return the Filter Plate and the Collection Plate to the manifold.
- Apply vacuum at -12 inches Hg for an additional 10 min.

- Add Wash Buffer.
- Centrifuge at 5,600-6,000 x g for 10 min
- Stand the Filter plate on a clean paper towel at room temperature for 5 min.

STEP 7. Plasmid Elution:

- Add Elution Buffer or ddH₂O to the Filter Plate. Stand for 3 min.
- Close the manifold valve. Turn on the vacuum source to build up a vacuum to -12 inches Hg.
- Open the manifold valve to apply vacuum to elute plasmid.
- Alternative:** If the consistent volume of elutes are recommend, use centrifuge protocol to process this elution step. (STEP 7-A-7-D).

- Add Elution Buffer or ddH₂O to the Filter Plate. Stand for 3 min.
- Centrifuge to elute plasmid.

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